


G-Series



711 Hulman St.
PO BOX 2128
TERRE HAUTE, IN 47802
812-235-6167
FAX 812-234-6975

DUCTLESS FUMEHOOD INSTALLATION AND OPERATION

- PARTS LIST:**
- Combined: Blower Unit Housing (A) and Filter Housing (B)
 - AC Line cord (C) (packaged inside Filter Housing) (B)
 - Enclosure Assembly (D)
 - Filters (Pre-filter & Carbon or Hepa) (E)

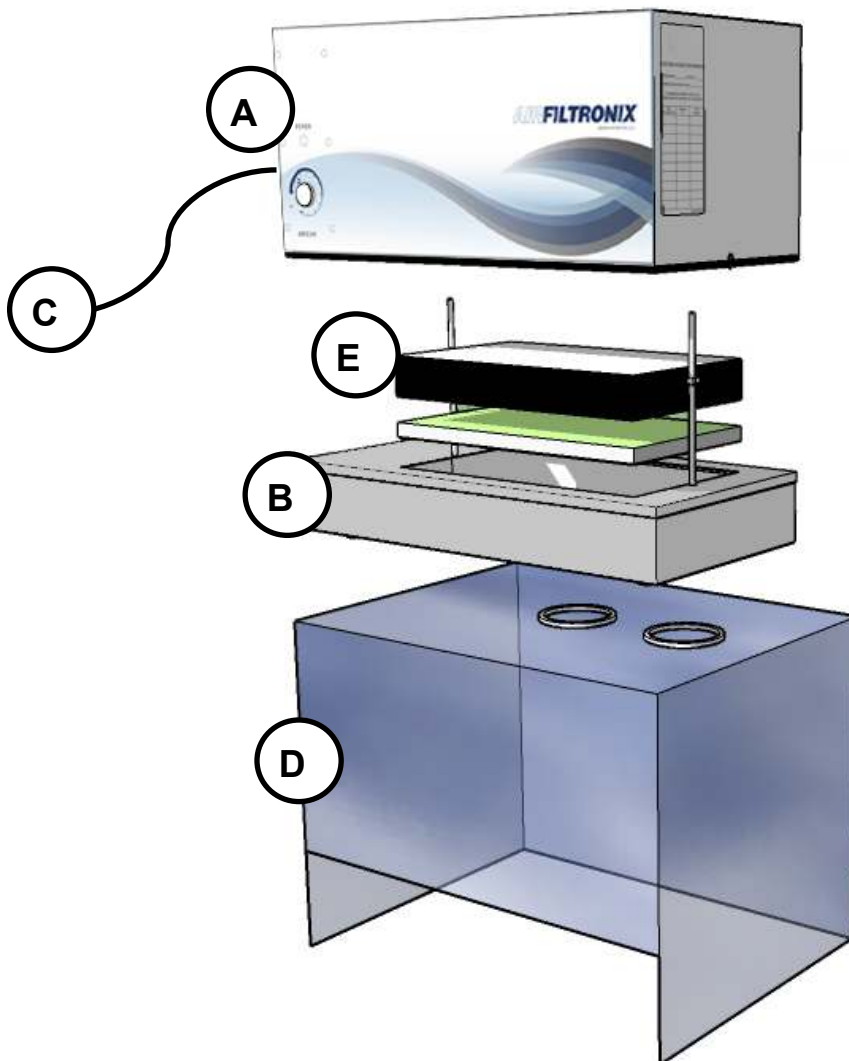

Remove section A from B by
unsnapping the 2 latches.

Install Filters: Pre-Filter first, then Car-
bon and/or Hepa with gasket down.

After filters are secured with bracket,
attach section A to B and place blower
unit on enclosure.

Filter Installation Guidelines: The Blower / Filter housing is designed to accommodate a variety of filter types which should be arranged accordingly: Pre-filters should be installed first, gasket side down. Activated carbon filters or other chemical absorbing or Hepa filters should be installed next, gasket side down. The filter installation / replacement history should be maintained using the card supplied in the clear pouch on the side of the blower. Tighten down both sides of the compression bar evenly with the wingnuts until the bar bends upward approximately 1/4 inch in the middle.



Once the unit is plugged into a properly grounded receptacle, turn the knob clockwise to activate the blower.

Safety Procedures: Turn off the blower and remove the line cord before replacing filters or cleaning the system. Wear OSHA approved gloves, safety glasses, and filtered breathing mask during maintenance procedures. Used filters should be placed in appropriate containers for disposal in accordance with local, state, and federal regulations.

Maintenance – Enclosure Cleaning: There are no user serviceable parts inside the blower unit housing. The filters should be checked for system performance and replaced on a regular basis. The surfaces of the enclosure can be cleaned with a solution of mild soapy water using a damp cloth or Windex™ type glass cleaner.

Limited Warranty:

See the Glas-col warranty policy located under the Warranty section on the Glas-Col website at www.glascol.com

Returns:

Call or fax Customer Service for a Return Goods Authorization (RGA) number before returning a Glas-Col product. Reference the RGA number on the shipping box and on a written description of the problem. A 20% restock charge of the net price is charged for all standard product returned to stock.

Glas-Col, LLC

711 Hulman Street, P.O. Box 2128, Terre Haute, IN 47802-0128

Phone. (812) 235-6167, Fax. (812) 234-6975

E-mail: pinnacle@glascol.com

Web: www.glascol.com