



HOTSPOT Heat Lamp

108A HL250-12



User Instructions

The HOTSPOT Heat Lamp assembly is shipped with the bulb removed. Screw the bulb into lamp socket and install the guard. Place three of the wire arms inside the rim of the shade then press fourth arm into position on the rim. (Note: for safety and comfort, guard must be attached whenever lamp is used).

To remove and Replace the Bulb, Unplug the power cord. Press any arm of the guard inward and pull down to remove guard from shade. Unscrew and replace the bulb - use 250W heat lamp only. Replace the guard.

To operate the device, plug the cord from the heat lamp into the provided variable control and then plug the control into a 120V outlet. Turn on the switch and adjust the knob to vary the intensity of the bulb. If the bulb does not light, rotate the switch on top of the heat lamp until the bulb illuminates and leave in the ON position. Caution: The housing of a heat lamp gets hot after a period of use. Use care when redirecting the position of the lamp when it is ON, or has just been turned OFF.

Spare Parts: 250-Watt infrared lamp (R64 574)

Limited Warranty:

See the current Glas-Col warranty policy located under the General Sales Policy on the Glas-Col website at www.glascol.com

LIMITATION OF WARRANTIES

APART FROM SUCH WRITTEN STATEMENT OF WARRANTY, THERE ARE NO WARRANTIES, EXPRESSED, IMPLIED OR STATUTORY, WHICH EXTEND BEYOND THE DESCRIPTION OF PRODUCTS ON THE FACE HEREOF.

Glas-Col products are intended only for legal and legitimate purposes in commercial laboratory settings.

Glas-Col reserves the right to make product refinements without prior notice.

Glas-Col, LLC

711 Hulman Street P.O. Box 2128, Terre Haute, IN 47802 0128 USA

Phone: (812) 235-6167 • Fax: (812) 234-6975

Email: pinnacle@glascol.com

www.glascol.com